

## (Re)starting the Mythen IOC

1. Log into mcspc01 as "rixs". This is the computer connected to the monitor on the upper left.
2. 'restartioc' to (re)start the IOC
3. 'restartmcs' to (re)start the MCS
4. 'start\_epics\_27mcs' to bring up the MEDM window for the MCS
5. cd /net/s27dserv/xorApps/epics/synApps\_5\_5/ioc/30idbMCS/30mcsApp/op/mythen/bin
6. ./start\_mmi\_dectris
7. Back in spec... 'mythen\_on'